


GXK-012

動物皿

取扱説明書

適用機種

- GX-K シリーズ全機種
- GF-K シリーズ全機種
- GX-M シリーズ全機種
- GF-M シリーズ全機種

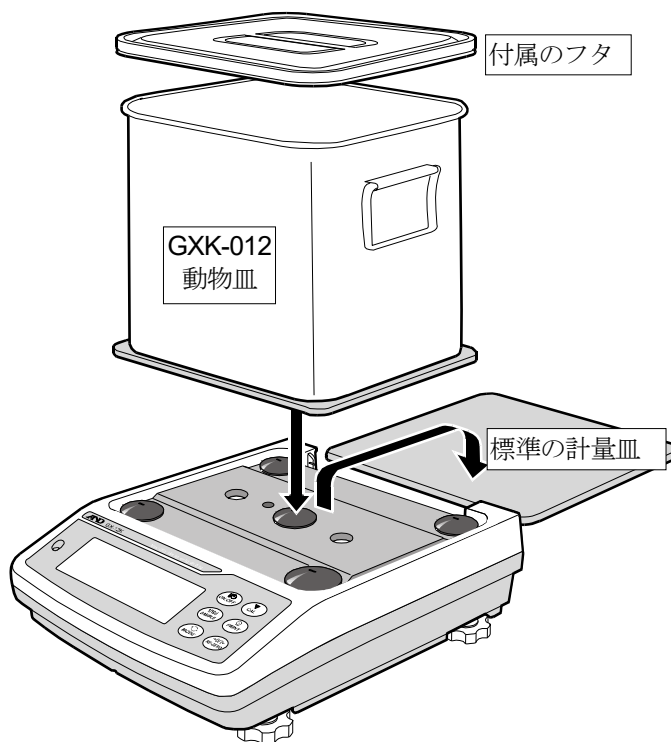

使用方法

準備

- 標準の計量皿を外し、GXK-012 動物皿に交換します。

計量

- RE-ZERO** キーを押し、表示をゼロにします。
- 動物を入れ、計量します。

注意

- 天びんのホールド機能（動物計量）を使用すると計量値の平均値を固定表示できます。
- 動物皿を使用した時はひょう量から約 1.5 kg 引いた値まで測定できます。
- 付属のフタは必要な場合に使用してください。
- 感度調整をする時は標準の計量皿を利用してください。
- 動物皿を載せた状態で感度調整し、その後標準の計量皿を載せると **-E** 表示になります。標準の計量皿に交換する場合は標準の計量皿を載せた状態で再度感度調整してください。


GXK-012

Animal Weighing Bowl

Instruction Manual

Models

- GX-K series
- GF-K series
- GX-M series
- GF-M series

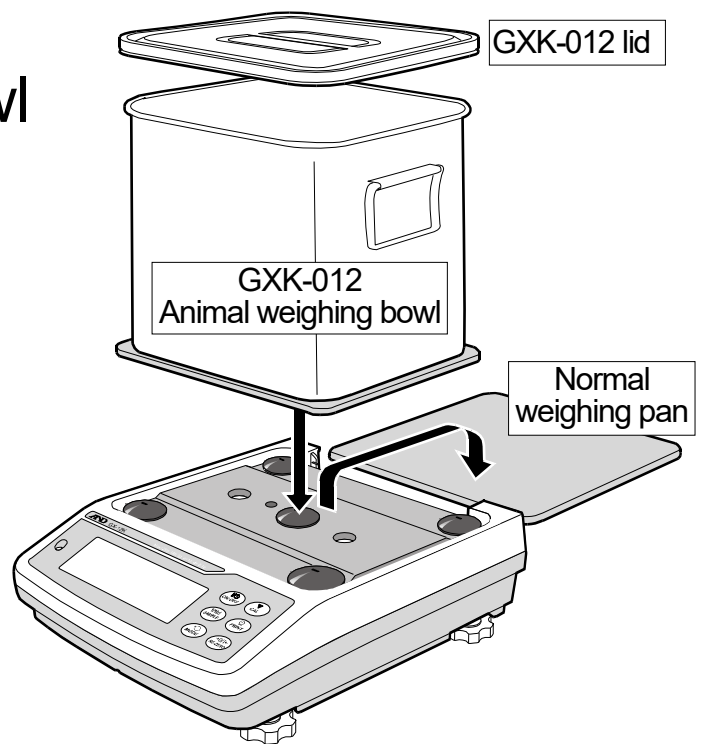

Weighing

Setup

1. Replace the normal weighing pan with the animal weighing bowl (GXK-012).

Measurement

2. Press the **RE-ZERO** key to cancel the weight (tare).
3. Place a sample animal in the bowl to weigh it.

Caution

- When using the animal weighing mode (Hold function), the average weight of the animal can be displayed.
- When using this option, the weighing range that can be used is an approximately 1.5 kg less than the weighing capacity.
- If necessary, use the GXK-012 lid.
- When performing sensitivity adjustment with the balance, use the normal weighing pan.
- If performing sensitivity adjustment with the balance that has the animal weighing bowl on it and then replacing it with the normal weighing pan, **-E** is displayed. When using the normal weighing pan again, re-perform sensitivity adjustment with the balance that has the normal weighing pan on it.

A&D

A&D Company, Ltd.