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Plastic Storage Container For SV Series Vibro Viscometer

AX-SV-39

Instruction Manual

1. Introduction

The AX-SV-39, a plastic storage container, is used to store sample fluids for an SV series Vibro Viscometer. It was designed to have an excellent sealing ability and should not leak if handled normally.

The viscosity of the sample stored in the container can be measured as it is.


Composition:

Plastic storage container (120 mL max.)	20 pcs
Material	
Container / lid	PP (Polypropylene)
	Operating temperature: 120°C max.
Inner lid	LDPE (Low density polyethylene)
	Operating temperature: 80°C max.

2. When used for measurement


An appropriate sample amount for measurement is approximately 100 mL.

(The volume up to a position 47 mm above the bottom of the plastic storage container corresponds to 100 mL.)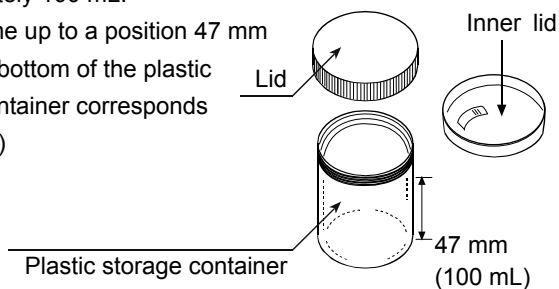


3. How to use

■ Place the plastic storage container on the table.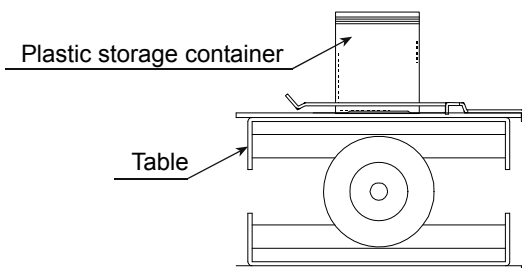

■ When the plastic storage container is to be used, be sure to raise or remove the protector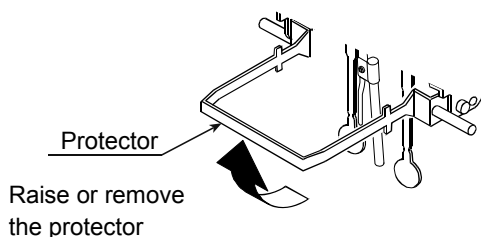


■ Use much care so that the sensor plates do not come into contact with the inner wall of the plastic storage container. If they are in contact, a stabilized display of the viscosity value is not obtained and the viscosity can not be measured.

■ Adjust the position of the plastic storage container so that the distance between the inner wall of the container and the right sensor plate is the same with the distance between the inner wall and the left sensor plate. If either one is closer to the wall, a measurement error will occur.

■ Adjust the table height so that the center of the narrow part of the sensor plates is at the sample surface.

At this time, use the surface locator plate as a guide. The surface locator plate has been secured in position so that the tip of the surface locator plate comes into contact with the sample surface. The position of the surface locator plate can be changed. (Refer to "12. TROUBLESHOOTING - How to remove and attach the surface locator plate" of the viscometer instruction manual.)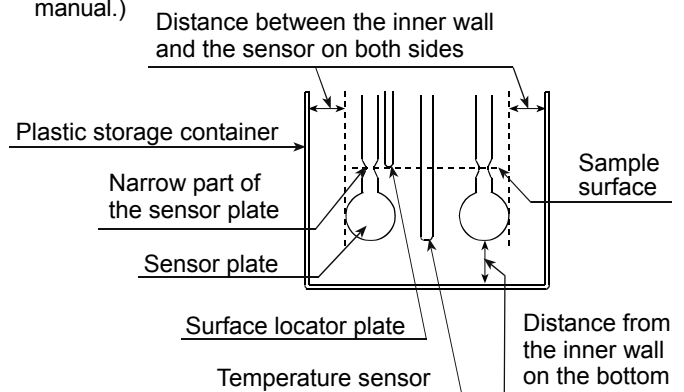

4. Measuring the viscosity

The viscometer has been calibrated using the accessory sample cup (35 mL) when shipped. The distance between the inner wall and the sensor plates when the accessory sample cup is used, differs from the distance when the plastic storage container is used. This causes a difference in the sensor plate's detection capability, thus causing a difference in the viscosity measured.

Therefore, to obtain an accurate viscosity value using the plastic storage container, it is recommended that calibration be performed using a fluid with a known viscosity value which is close to the sample viscosity. (Refer to "6. VISCOSITY CALIBRATION" of the viscometer instruction manual.)

5. Maintenance

Clean the plastic storage container as necessary. If cleaning is not sufficient, a measurement error may occur due to contamination.