


FXi-12

動物計量皿

取扱説明書

適用機種

FX-1200i / 2000i / 3000i

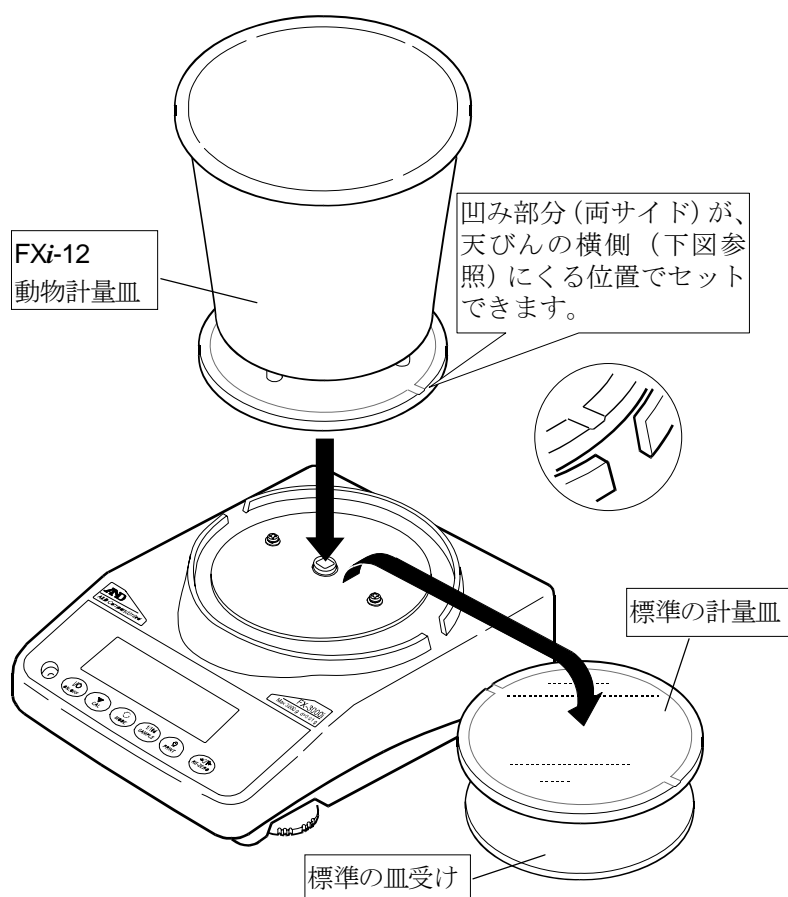

使用方法

準備

- 1 標準の計量皿と皿受けを外し、FXi-12 動物計量皿に交換します。

計量


- 2 **RE-ZERO** キーを押し、表示をゼロにします。
- 3 動物を入れ、計量します。

注意

- 天びんのホールド機能(動物計量)を使用すると計量値の平均値を固定表示できます。詳細は、天びん本体の取扱説明書を参照してください。(内部設定 *bRSFnc Hold*、「環境・表示の解説」を参照願います。)
- 動物計量皿を使用した時はひょう量から約200g引いた値まで測定できます。
- キャリブレーションをする時は標準の計量皿を利用してください。
- 動物計量皿を載せた状態でキャリブレーションし、その後標準の計量皿を載せると **-E** 表示になります。標準の計量皿に交換する場合は標準の計量皿を載せた状態で再度キャリブレーションしてください。
- FX-120i / 200i / 300i には、動物計量皿(FXi-12)は使用できません。(動物計量皿の自重(約200g)により、天びんの使用可能なひょう量が無くなります。)

FXi-12

Animal Weighing Bowl Instruction Manual

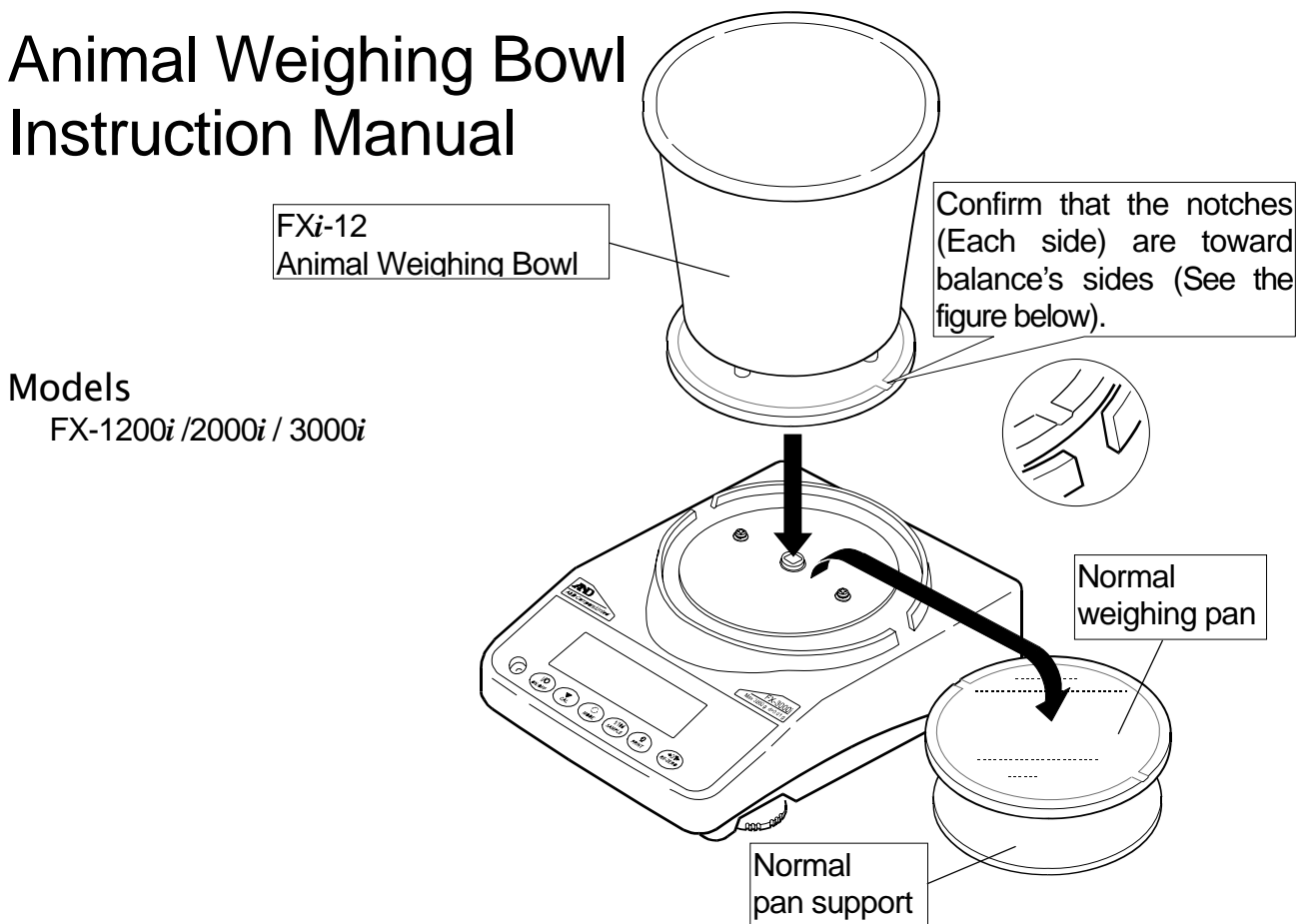

Weighing

Setup

- 1 Replace the normal weighing pan and pan support with the animal weighing bowl (FXi-12).

Measurement

- 2 Press the **RE-ZERO** key to cancel the weight (tare).
- 3 Place a sample animal in the bowl to weigh it.

Caution

- When using the animal weighing mode (Hold function), the average weight of the animal can be displayed. For details, refer to the instruction manual of the balance. (Refer to the function *bASFnc Hold*, Description of the Class “Environment, Display”)
- When using this option, the weighing range that can be used is an approximately 200 g less than the weighing capacity.
- When calibrating the balance, use the normal weighing pan.
- When calibrating the balance with the animal weighing bowl and replacing with the normal weighing pan, **-E** is displayed. If the normal weighing pan is used, re-calibrate the balance with it.
- The animal weighing bowl can not used with FX-120i / 200i / 300i. (The balance's weighing capacity that is available is used by the animal weighing bowl's weight (Approx. 200 g).)