

FXi-12

動物計量皿

取扱説明書


適用機種

FX-1200i / 2000i / 3000i / 5000i
FZ-1200i / 2000i / 3000i / 5000i

FX-1200i WP / 2000i WP / 3000i WP
FZ-1200i WP / 2000i WP / 3000i WP

FX-1202 / 2202 / 3202 / 5202
FZ-1202 / 2202 / 3202 / 5202

FX-1202WP / 2202WP / 3202WP
FZ-1202WP / 2202WP / 3202WP

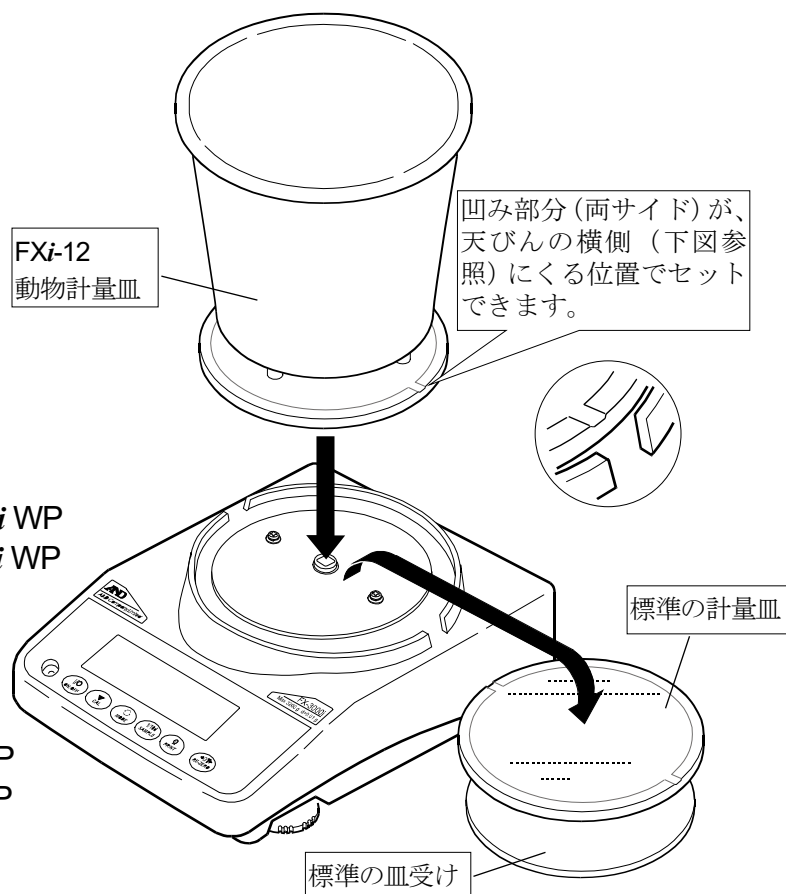

使用方法

準備

- 1 標準の計量皿と皿受けを外し、FXi-12 動物計量皿に交換します。

計量

- 2 **RE-ZERO** キーを押し、表示をゼロにします。
- 3 動物を入れ、計量します。

注意

- 天びんのホールド機能(動物計量)を使用すると計量値の平均値を固定表示できます。詳細は、天びん本体の取扱説明書を参照してください。(内部設定 *bR5Fnc Hold*、「環境・表示の解説」を参照願います。)
- 動物計量皿を使用した時はひょう量から約200g引いた値まで測定できます。
- 感度調整をする時は標準の計量皿を利用してください。
- 動物計量皿を載せた状態で感度調整し、その後標準の計量皿を載せると **-E** 表示になります。
標準の計量皿に交換する場合は標準の計量皿を載せた状態で再度、感度調整してください。
- FX-i / FZ-i、FX-i WP / FZ-i WP、FX / FZ、FX-WP / FZ-WP の最小表示が1mgの機種には、動物計量皿(FXi-12)は使用できません。(動物計量皿の自重(約200g)により、天びんの使用可能なひょう量がなくなります。)

FXi-12

Animal Weighing Bowl Instruction Manual

FXi-12
Animal Weighing Bowl

Confirm that the notches
(Each side) are toward
balance's sides (See the
figure below).


Models

FX-1200i / 2000i / 3000i / 5000i
FZ-1200i / 2000i / 3000i / 5000i

FX-1200i WP / 2000i WP / 3000i WP
FZ-1200i WP / 2000i WP / 3000i WP

FX-1202 / 2202 / 3202 / 5202
FZ-1202 / 2202 / 3202 / 5202

FX-1202WP / 2202WP / 3202WP
FZ-1202WP / 2202WP / 3202WP

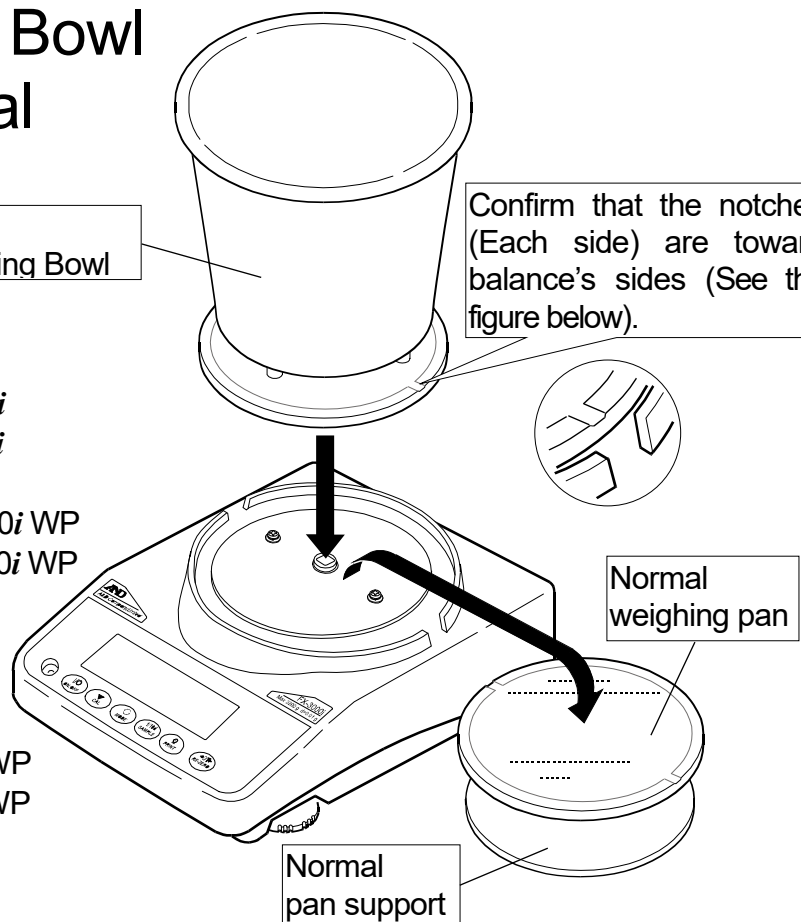

Weighing

Setup

- 1 Replace the normal weighing pan and pan support with the animal weighing bowl (FXi-12).

Measurement

- 2 Press the **RE-ZERO** key to cancel the weight (tare).
- 3 Place a sample animal in the bowl to weigh it.

Caution

- When using the animal weighing mode (Hold function), the average weight of the animal can be displayed. For details, refer to the instruction manual of the balance. (Refer to the function *bASFnc Hold*, Description of the Class “Environment, Display”)
- When using this option, the weighing range that can be used is an approximately 200 g less than the weighing capacity.
- When performing sensitivity adjustment with the balance, use the normal weighing pan.
- If performing sensitivity adjustment with the balance that has the animal weighing bowl on it and then replacing it with the normal weighing pan, **-E** is displayed. When using the normal weighing pan again, re-perform sensitivity adjustment with the balance that has the normal weighing pan on it.
- The animal weighing bowl can not used with FX-i / FZ-i, FX-i WP / FZ-i WP, FX / FZ, FX-WP / FZ-WP models with a readability of 1 mg. (The balance’s weighing capacity that is available is used by the animal weighing bowl’s weight (Approx. 200 g).)