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
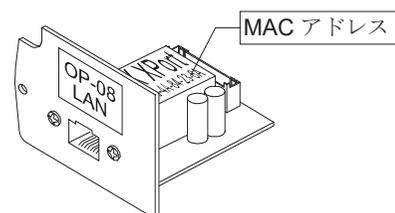
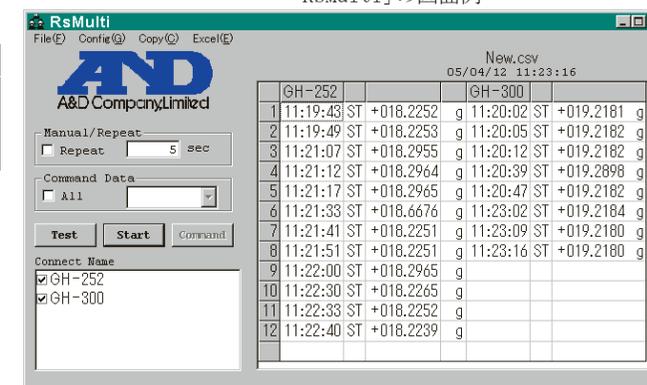
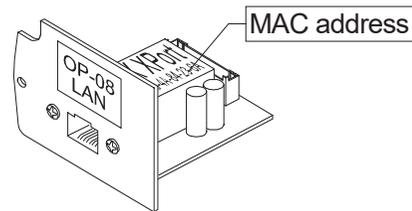
Unpacking the GH-08

GH-08 board
Ethernet Interface

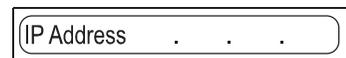AC adapter labels
Affix them on the AC adapter.

IP address label (5 labels)
Write a preset IP address and affix it on the balance.

Protective cover (5 labels)
Affix it on the IP address label.

Installing the GH-08

Step 1 Record the MAC address.
The IP address can be specified with it.

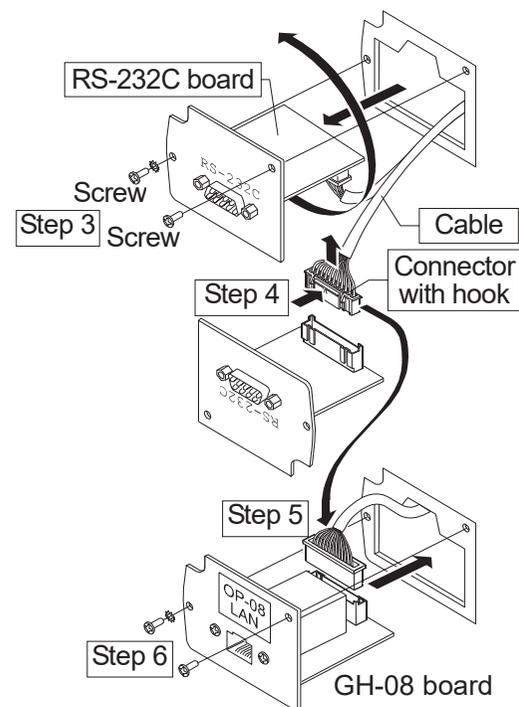
Step 2 Remove the AC adapter from the balance.

Step 3 Remove two screws and pull out the RS-232C board gently. Do not pull the cable excessively.

Step 4 Push the hook aside and remove the connector from the RS-232C board.

Step 5 Insert the connector into the GH-08 board.

Step 6 Fix the GH-08 board with two screws.

Wiring to the Computer

Caution

- Do not use the GH-08 in an important network.
A&D and dealers make no warranties if the GH-08 creates a network error.
- Contact the network manager before connecting the GH-08 to a network.
- Preset the IP address and subnet mask to the GH-08 one by one.
The same IP address 172.16.100.2 is stored at the factory. Do not duplicate an IP address.
- The IP address can not be returned to factory setting.
We recommend that you write the IP address on the supplied label.

Preparation

Please download the following software from A&D website:

<https://www.aandd.jp/products/software/software.html>

- Setting procedure "WinCT-Plus Instruction Manual"
- IP address setting software "Device Installer"
- Windows communication software "WinCT-Plus"

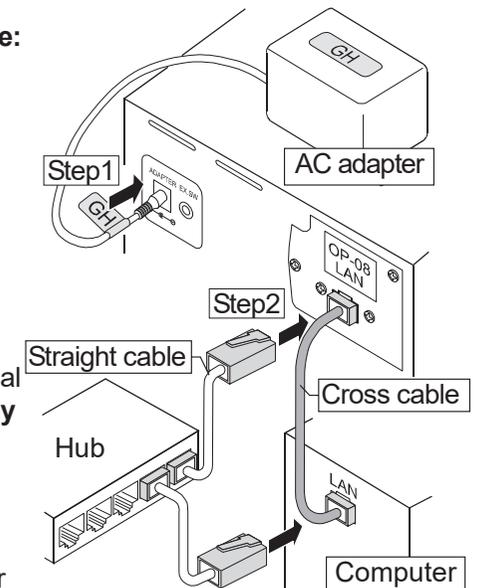
Step 1 Connect the AC adapter supplied with the balance to the balance.

Step 2 Connect the GH-08 and the computer with a cross cable directly or connect them with hub and straight cables.

Step 3 Input the IP addresses and a subnet mask of the GH-08 and computer. Refer to the WinCT-Plus Instruction Manual for the setting procedures. **Communication errors may occur with the wrong operation.**

Step 4 Write the IP address on "the IP address label" and affix it on the balance. You can see the address easily.

Step 5 Install the data acquisition software "RsMulti" into the computer. Refer to the WinCT-Plus Instruction Manual for the installation procedure.

The Outline of the Data Acquisition Software


Run the data acquisition software "RsMulti". Menu: [Start] → [Program] → [A&D WinCT-Plus] → [RsMulti]. Refer to the [A&D WinCT-Plus] → [Manual] for the operation.

- The software can acquire data from multiple devices connected with a LAN or RS-232C.
- The software can control these devices with commands.
- The software can acquire data transmitted from devices.
Example: When pressing the **PRINT** key of the balance, data is output and is acquired with the computer.
- The stored data can be used with EXCEL. (Install the Microsoft EXCEL before use.)

Example

Two balances are connected.

	IP address	Subnet mask
Computer	172.16.100.1	255.255.0.0
GH-252	172.16.100.2	
GH-300	172.16.100.3	

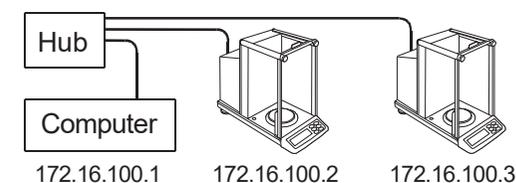A sample window of the "RsMulti"