AD-5701BB

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1WMPD4005680

© 2025 A&D Company, Limited.

No part of this publication may be reproduced, transmitted, or translated into any language in any form by any means without the written permission of A&D Company, Limited.

The contents of this manual and the specifications of the instrument covered by this manual are subject to change for improvement without notice.

Other trademarks and trade names are those of their respective owners.

Introduction

This manual describes how the product, the AD-5701BB (3-Channel 100-Hour Digital Timer), works and how to get the most out of the product. Read this manual thoroughly before using the product and keep the manual to hand for future reference.

Safety Precautions

All safety messages are in accordance with the ANSI Z535.4 (American National Standard Institute: Product Safety Signs and Labels) standard. **WARNING DEFINITIONS**

The warnings described in this manual mean the following:

	This symbol indicates a potentially hazardous situation that, if not avoided, could result in personal injury or property damage. It may also be used to alert the user about unsafe practices.
--	---

When using the AD-5701BB, the following safety precautions should always be followed.

Caution


Internal servicing or adjustment of the product must be performed by a qualified person.

Precautions for Use

- This product is a precision instrument and should be handled with care. Please be careful not to drop the product or subject it to strong impacts, as this may cause damage or malfunction.
- When attaching a magnet to a metal surface, there is a risk of scratching the metal surface. Therefore, please test on an inconspicuous area before use.
- Do not use the product in places where it will be exposed to direct sunlight for a long time, in a closed car, or near a heater such as a stove. The operating temperature range of this product is 0 to 40°C. Use of the product beyond this temperature range may cause the product to malfunction.
- Avoid using the product in places where there can be sudden changes in temperature, high temperature, high humidity, or high dust content.
- Avoid suddenly moving the product from a hot environment to a cold one or vice versa. Sudden temperature changes may cause water droplets to form inside the product.
- Do not place the product in a humid or dusty place, as this may cause the product to malfunction. The operating humidity range of this product is 10 to 90% RH. Water accidentally entering the product may cause a malfunction. Do not place objects filled with water around this product.
- Places with strong magnetic or electric fields (TVs, induction cookers, microwave ovens, etc.) may affect the device. Please avoid using the product in such places.
- To prevent hazards, do not use the product in the presence of flammable gases.
- The product is not waterproof. Do not use the product in places where it will


be exposed to rain or water.

Part Names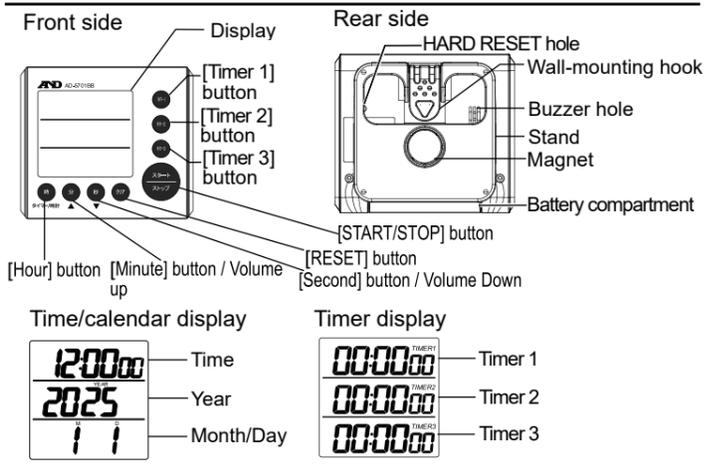

Placement

This product can be hung on a wall (screws not included), hung around the neck by putting a strap through the wall mounting hook (strap not included), attached to a metal surface with a magnet, placed upright on a tabletop using the stand, or clipped to a pocket or belt.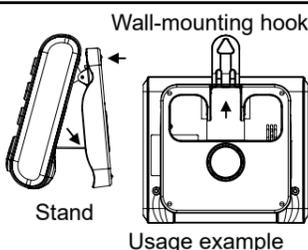


Installing / Replacing Battery

Please insert the battery following the correct procedure below. When the battery runs low, the display may become difficult to read, or the device may malfunction or become unresponsive. If this occurs, please replace the battery.

Since the battery compartment is located under the stand, you will need to lift the stand to access it.

Note that replacing the battery will erase all stored information such as the time, date, and timer settings, so these will need to be set again.

- Remove the battery cover on the back of the main unit.
- Remove the old battery.
- Install a new AAA battery, oriented correctly so that the polarities match those shown inside the compartment.
- Put the battery cover back on.

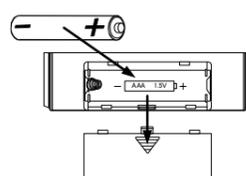Caution

- The included battery is for testing and its life may be short.
- If the product is not going to be used for a long time, remove the battery and store in a safe place.
- The battery life described in the specifications is based on when a new alkaline battery is used at an ambient temperature of 25 °C and the buzzer is used 3 times a day (approx. 90 seconds). Depending on the ambient temperature and the frequency of buzzer use, the battery life may become extremely short.
- Make sure to use an AAA (cell) battery only.
- Be sure to insert the battery with the polarities (+/-) in the correct orientation. If the batteries are inserted with the polarities in the wrong orientation, not only will the product not work properly, but it may also cause the product to malfunction.
- Do not charge, short-circuit, disassemble, or expose the battery to fire as this may cause the battery to explode or leak.
- Keep batteries out of the reach of children.

Operations (Time and Date Settings)

Switching between Time & Date Display and the Timer Display

This product can show either the Time & Date display or the Timer display. The Time & Date display shows the current time and date (year, month, and day). The Timer display shows three timers. To switch between the displays, hold down the [Hour] button for 2 seconds or longer.

Setting the Time and Date

- While the time and date are displayed, press the [Minute] and [Second] buttons

at the same time. The time display blinks and the device enters the Date setting mode.

- Press the [Hour] button to advance the time by 1 hour. Hold the button down to advance the time quickly.
 - Press the [Minute] button to advance the time by 1 minute. Hold the button down to advance the time more quickly.
 - Press the [Second] button to advance the time in 10-second increments. Hold the button down to advance the time more quickly.
- Press the [START / STOP] button once to change the year setting. The year display will start blinking.
 - Press the [Minute▲] button to advance the date by 1 year. Hold the button down to advance the date more quickly.
 - Press the [Second▼] button to set the date back by 1 year. Hold the button down to set the date back more quickly.
 - Press the [RESET] button to reset the year display to the default, 2025. The setting range is from 2025 to 2099.
 - Press the [START / STOP] button once again to change the month setting. The month display will start blinking.
 - Press the [▲] button to advance the date by 1 month. Hold the button to advance the date more quickly.
 - Press the [▼] button to set the date back by 1 month. Hold the button to set the date back more quickly.
 - Press the [RESET] button to reset the month display to the default, 1.
 - Press the [START / STOP] button once again to change the day setting. The day display starts blinking.
 - Press the [▲] button to advance the date by 1 day. Hold the button to advance the date more quickly.
 - Press the [▼] button to set the date back by 1 day. Hold the button to set the date back more quickly.
 - Press the [RESET] button to reset the day display to the default, 1.
 - Press the [Start / Stop] button once again to exit time and date setting and return to the Time & Date display.
 - Note: The display goes back to the Time & Date display if buttons are not touched for 2 minutes during time and date setting.

Switching between 12 and 24-Hour Formats


When the time and date are displayed, pressing the [Hour] button switches between the 12- and 24-hour formats. 'AM' or 'PM' is displayed in the 12-hour format.

Operations (Timer Settings)

Timer Setting

If the display shows the time and date, hold down the [Hour] button for 2 seconds or longer to switch it to the Timer display. A separate timer can be set for each of the three channels: Timer 1, Timer 2, and Timer 3. Each timer can be set to up to 99 hours 59 minutes and 59 seconds.

- When the Timer display is shown, hold down the [Timer] button of the channel you want to set for 2 seconds or longer. The time in that channel in the display starts blinking. The image on the right is an example for Timer 1. The [Hour] button advances the timer by 1 hour, the [Minute] button by 1 minute, and the [Second] button by 1 second. Press and hold each button to advance the time more quickly. Press the [RESET] button to reset the timer to 00:00:00.
- After setting the timer, press the [Timer] button of the channel you are setting. The display stops blinking and the setup is complete. SET INDENT TO MATCH (1), (2) If no buttons are pressed for 2 minutes or longer during the setup, the Timer 1 display stops blinking and the setup is stopped. To set up a timer again, restart from step 1 above.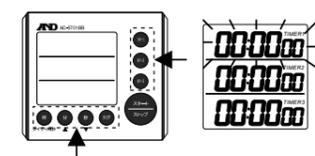

How to Use the Timer

The three channels (Timer 1, Timer 2, and Timer 3) can be used independently. Alternatively, all three channels can be started simultaneously.


o To Use Timer 1, Timer 2, and Timer 3 Independently

- When the Timer display is shown, set the timer for the channel you want to use by following the steps for 'Timer Setting.'
- Press the [Timer] button of the channel which the timer has been set to start countdown. Pressing the [Timer] button of a channel without a timer set starts count-up and displays a blinking 'TIME'S UP' message on the left side.
- Press the same [Timer] button as done in step 2 to stop countdown/count-up.

o To Start Timer 1, Timer 2, and Timer 3 Simultaneously

- When the Timer display is shown, follow the steps for 'Timer setting' to set a timer for Timer 1, Timer 2, and Timer 3.
- Press the [START / STOP] button to start the countdown of Timer 1, Timer 2, and Timer 3 simultaneously. A count-up starts for channels that do not have a timer set and a blinking 'TIME'S UP' is shown on the left side.
- Press the [START / STOP] button while the countdown/count-up is being performed to stop the countdown/count-up of Timer 1, Timer 2, and Timer 3. Press the [Timer] button of a channel while its countdown/count-up is being performed to stop the countdown/count-up of that channel only.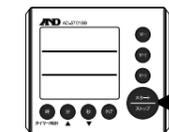

Operation When the Set Time of the Timer Is Reached

- When the set time is reached, 'TIME'S UP' is displayed on the left side of the channel that has finished. Then, the count-up starts and the buzzer beeps for approximately 30 seconds. The buzzer types vary depending on each channel as follows.
 - Timer 1: beep, beep, beep...
 - Timer 2: beep-beep, beep-beep, beep-beep...
 - Timer 3: beep-beep-beep, beep-beep-beep, beep-beep-beep...
- * The buzzer automatically stops after approximately 30 seconds but, the count-up does not stop.
- * A buzzer stops beeping when the buzzer for another channel starts beeping.
- To stop a buzzer, press any button ([Hour]/[Minute]/[Second]/[RESET]) other than the [START / STOP] button. The buzzer for the channel that is beeping stops, and the count-up stops as well. Other channels continue their countdown/count-up. Please note if you press the [Start/Stop] button while the buzzer is beeping, the timers for all three channels will stop along with the buzzer.

Using the Same Timer Settings Again (Reload)

- Press the [RESET] button to display '00:00:00' for channel(s) for which the timer operation has stopped.
- Press the [Timer] button of a channel which displayed '00:00:00' in step 1 to see the previous setting value. If all channels displayed '00:00:00', press the [Start/Stop] button to see the previous setting values for all channels.
- Press the [Timer] button of the applicable channel again to start the countdown.

Volume Adjustment

While the timer is displayed, pressing the [Minute▲] button increases the volume and pressing the [Second▼] button decreases it. If no operation is performed for 3 seconds, the display will return to the timer screen.

Hard Reset

If this product stops working properly, it can be reset by performing the following two steps.

- Remove the battery. Wait for 30 seconds and reinsert the battery
 - Insert something pointed (e.g. paperclip) into the HARD RESET hole on the back of the product
- SET INDENT TO MATCH (1), (2) When a hard reset is done, the time, date, and timer setting values are all reset to the factory settings. Please set them again.

Maintenance

To clean the product, first dampen a soft cloth and then wipe gently. Do not use thinner, benzene-like volatile solvents, or abrasives.

Troubleshooting

Nothing is displayed	Make sure there is a battery in the battery compartment. If there is, it may have been depleted. Replace it with a new battery.
The display is too dim	The remaining battery life is short. Replace the battery with a new one.
The product does not work properly	The internal circuitry may have stopped working. Hard reset the timer. When the battery is low, the product may not work properly, so please replace the battery.

Specifications

Display style	: Switching between 'Time & Date (year, month, day)' and 'Three-channel Timer'
Alarm setting range	: 00:00:00 to 99:59:59
Number of channels	: 3
Timer functions	: Count-up, Count-down
Buzzer sound duration	: 30 seconds
Buzzer volume	: Off / Lo : 60 dB at 10 cm / Hi : 75 dB at 10 cm
Time display	: Current time (hh:mm:ss), switches between 12-/24-hour formats, Current date (year/month/day)
Accuracy	: ±1 second/day
Battery	: 1 AAA (cell) battery
Battery life	: Approx. 1 year (when an alkaline battery is used and the alarm is operated for 90 seconds a day)
Dimensions	: 80 (W) x 70 (H) x 30 (D) mm
Weight	: Approx. 85g (incl. battery)
Operation environment	: 0 to 40°C, 10 to 90%RH (without condensation)
Storage environment	: -10 to 50°C, 10 to 95%RH (without condensation)
Items included	: Battery (for monitoring), instruction manual
Material	: Upper cover, lower cover, side panel, stand, battery cover, wall-mounting hook: ABS Buttons: Rubber Metal parts for the stand: SUS304 Screws: SUS430